# Scout<sup>®</sup> SKX Series Balances

## User Guide

EN

ES

FR

DE

IT

PT

HU

SV

DA

PL

CS

NL

# 1. SAFETY INFORMATION

## Definition of Signal Warnings and Symbols

- WARNING** For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
- CAUTION** For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
- Attention** For important information about the product.
- Note** For useful information about the product.

## Warning Symbols

General hazard

Electrical shock

## Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

## Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.


# 2. INSTALLATION

## 2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.


## 2.2 Releasing the Transportation Lock

On top of the balance, turn the pointer 90 degrees counter-clockwise to unlock.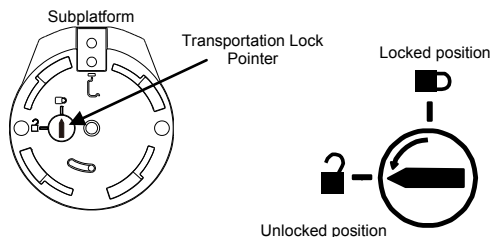

### 2.3 Installing the Weighing Pan

Balances with a rectangular pan are placed into the sub-pan as shown and rotated counter-clockwise until it locks. Round pans are placed straight down on sub-pan.

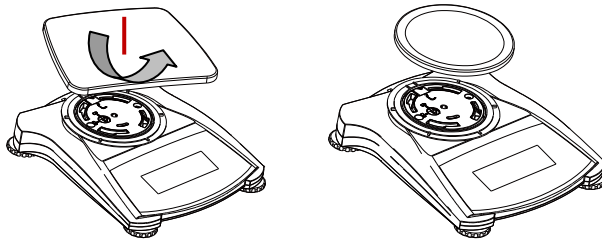
### 2.4 Connecting Power

#### AC Adapter Installation

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

**Attention:** Only use a AC adapter specified by OHAUS.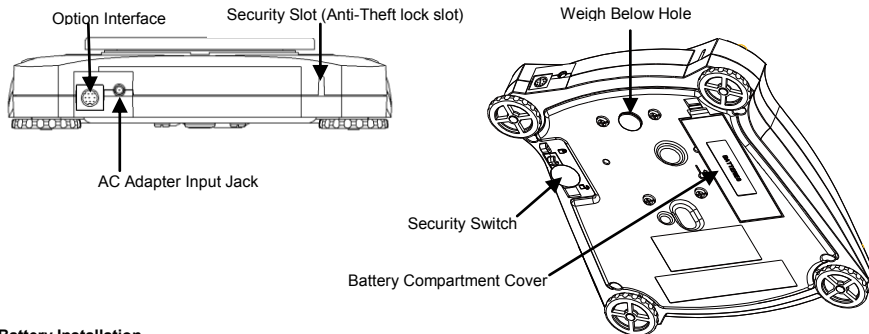

#### Battery Installation

Install the four "AA" batteries with polarity as shown in the battery compartment.

### 2.5 Initial Calibration


For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

## 3. OPERATION

### 3.1 Overview of Display, Home Screen

#### Button

#### Functions:

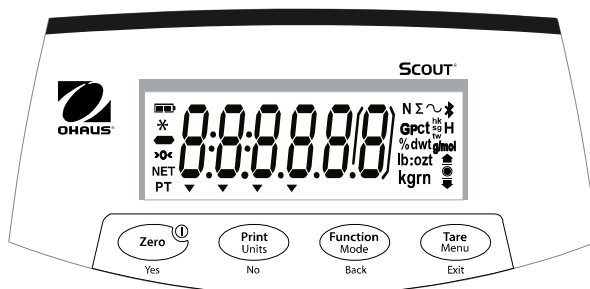Button	Zero Yes	Print Units No	Function Mode Back	Tare Menu Exit
Short Press	<b>Zero/On</b> Turns the balance on; If balance is On, sets Zero	<b>Print</b> Initiates print	<b>Function</b> Initiates an application mode	<b>Tare</b> Tare
Long Press	<b>Off</b> Turns the balance Off	<b>Units</b> Changes the weighing unit	<b>Mode</b> Changes the application mode	<b>Menu</b> Enters the User menu
Short Press (Menu Function)	<b>Yes</b> Accepts the current setting	<b>No</b> Advances to the next menu or setting	<b>Back</b> Moves to previous menu item	<b>Exit</b> Exits the User menu

### 3.2 Menu Structure

Cal	Setup	Mode	Unit	End
Span	Reset	Reset	Reset	
Lin	Filter	Weighing	g	
	AZT	Density	kg	
	StableRange	Mole Weighing	oz	
	Backlight		lb:oz	
	Auto Tare		lb	
	Auto Off		N	
End	End	End	End	

### 3.3 Using the Balance

Note: Before using any application, be sure the balance has been leveled and calibrated.

#### Weighing Application

1. If required, place an empty container on the pan and press **Tare**.
2. Add sample to the pan or container. The display shows the weight of the sample.

## 4. MAINTENANCE

### 4.1 Cleaning

**WARNING:** Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

**Attention:** Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

## EN-4

### 4.2 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website [www.ohaus.com](http://www.ohaus.com) to locate the Ohaus office nearest you.

## 5. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Specific temperature range: +10 °C to +40 °C
- Operating temperature range: +5 °C to +40 °C
- Relative humidity: 10% to 80% at 31 °C, decreasing linearly to 50% at 40 °C, non-condensing
- Mains supply voltage fluctuations: up to ±10% of the nominal voltage
- Installation category II
- Pollution degree: 2

### SPECIFICATIONS:

Model	SKX123	SKX222	SKX422	SKX622	SKX1202	SKX2202	SKX421	SKX621	SKX2201	SKX6201	SKX8200
Capacity (g)	120	220	420	620	1200	2200	420	620	2200	6200	8200
Readability (g)	0.001	0.01	0.01	0.01	0.01	0.01	0.1	0.1	0.1	0.1	1
Repeatability (Std. Dev.) (g)	0.002	0.01	0.01	0.01	0.02	0.02	0.1	0.1	0.1	0.1	1
Linearity (g)	0.003	0.01	0.01	0.02	0.03	0.03	0.1	0.1	0.1	0.2	1
Power	Standard AC Power Adapter – 5VDC 1A output, or 4 AA dry battery (not included)										

## 6. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> .
	AS/NZS 61000.6.1, AS/NZS 61000.6.3

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P/N 30269066 B © 2015 Ohaus Corporation, all rights reserved.
Ohaus Corporation